



Labels 100 x 100

⚠ WARNING

Contents under pressure.

Depressurize tank before removing eyewash assembly.

👁️ + PORTABLE EYEWASH/ DRENCH HOSE FILLING INSTRUCTIONS

1. Depressurize tank by activating the pressure relief valve on the out system.
2. Remove the eyewash assembly from the tank by removing the clamp.
3. Fill the tank with 42.5l (11.2 US gal) of clean water only (ISO 9001:2015 clause 7.1.4).
4. Add 100ml (3.38 US fl. oz) of disinfectant to the tank (ISO 9001:2015 clause 7.1.4).
5. Problem removal is essential, in order to keep an optimal safety status.
6. Pressure on the tank to 6.5 bar (95 PSI) - MAXIMUM - by way of the standard pressure relief valve (1.6 MPa or 23200 PSI) on the tank.
7. Turn the unit momentarily to purge air and any contaminants from systems.

Note: Water should be changed and replaced with fresh water and one cubic of bacteriostatic additive at least once every six months.

CAUTION: KEEP UNIT AWAY FROM HEAT

BROEN-LAB



						Description: Broen-Lab Portable Eyeshower				
						Media: Water				
						Comments: Portable Eyewasher, Vessel				
						Drawing type: Customer drawing				
Rev.	Pos.	Date	TM No.	Design by	Drawn by:	Change:	Designed by:	Drawn by:	Released on TM No.	Date (DD-MM-YY):
							MJO	MJO		18-06-25
BROEN-LAB						Format: A3		Drawing No.:		2737001100001
						This drawing is our property and cannot be given to a third party, copied or otherwise used without our permission. BROEN-LAB A/S				