


## Assembly, Operation and Maintenance Guide

2737001100001 / 2737002100001

### Assembly and Filling Instructions

- 1- **IMPORTANT:** Depressurize tank by activating pressure relief valve on the outlet stem assembly. Once unit is depressurized, activate flag handle on the eyewash to ensure unit is completely depressurized. Keep flag handle open. **Note:** Depressurize tank completely before proceeding to step 2.
- 2- Loosen and remove (2) bolts, nuts, clips, and the clamp from the tank, raise the eyewash assembly from the tank.
- 3- Fill tank with 37,5 L of clear clean potable water only. Then add entire bottle of bacteriostatic additive included with the unit. **DO NOT OVERFILL!**
- 4- Replace eyewash assembly ensuring that the eyewash assembly flag handle and the eyewash sign on the vessel are facing the same direction. Reclamp securely in place by tightening the nuts. Replace clips.
- 5- Using the Schrader air intake valve, pressurize unit to 6,5 bar. Be sure to use a clean air source. **DO NOT USE CO<sub>2</sub> OR ANY OTHER HIGH PRESSURE GAS! DO NOT CONNECT DIRECTLY TO ANY EXTERNAL AIR OR GAS SOURCE!**
- 6- Run unit momentarily to purge air and contaminants from system.

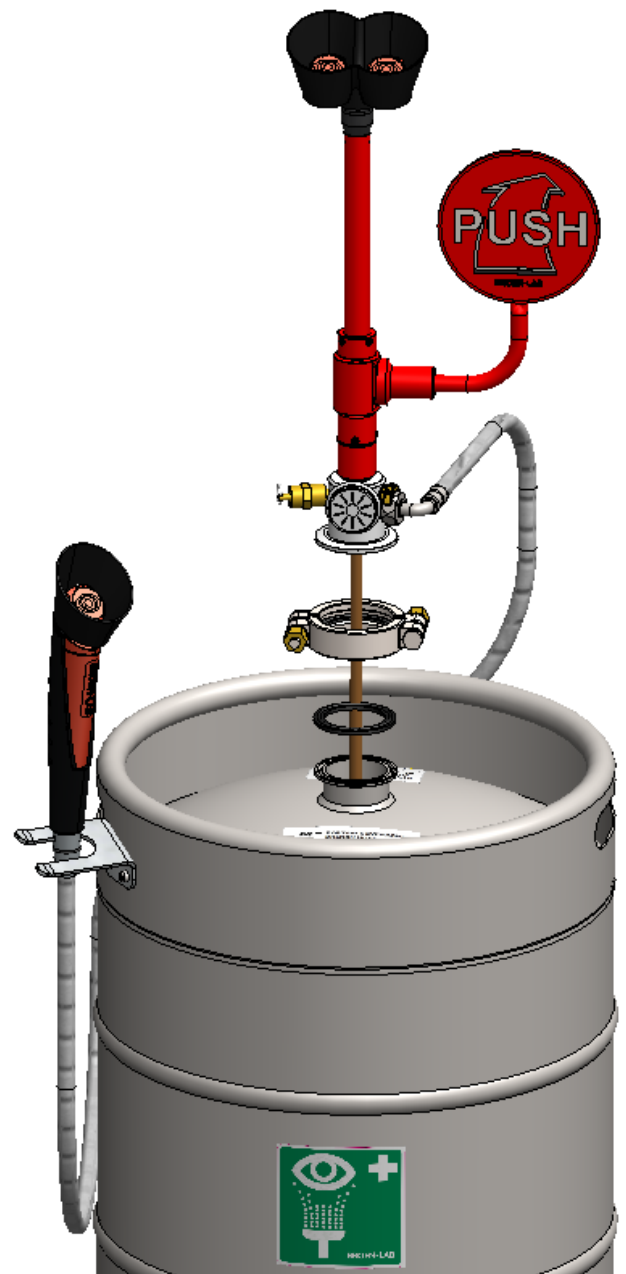

**Note:** Water should be drained and replaced with fresh water and one bottle of bacteriostatic additive at least once every six months.

**CAUTION: KEEP UNIT AWAY FROM HEAT!**

## ***Decommissioning of Unit***

- 1- **IMPORTANT:** Depressurize tank by activating pressure relief valve on the outlet stem assembly. Once unit is depressurized, activate flag handle on the eyewash assembly to ensure unit is completely depressurized. Keep flag handle open. **NOTE:** Depressurize tank completely before proceeding to step 2.
- 2- Loosen and remove (2) bolts, nuts, clips and the clamp from the tank. Raise the eyewash assembly from the tank.
- 3- Drain water from the vessel.
- 4- Replace eyewash assembly. Reclamp securely in place by hand tightening the nuts. Replace clips.

## ***General Notes***

The Pressurized transportable self-contained eyewash is designed to supply 2,5 L/min of water for 15 minutes in environments where a permanent water supply may not be available. Unit should be placed in a location where the spray nozzles are not less than 800 mm and no greater than 1200 mm above the surface on which the user stands. It shall also be positioned so that there are no obstructions on either side of the unit. Unit should be located on a surface which is flat, dry and out of direct sunlight.

USE OF THIS UNIT FOR ANY OTHER PURPOSE NOT INTENDED BY THE MANUFACTURER WILL VOID THE WARRANTY.

## ***Safety and Maintenance Instructions***

Read and follow all instructions for the eyewash station. Failure to do so may result in serious eye damage and vision loss.

Pressure relief valve should be checked to make sure it is working properly.

It is recommended that the eyewash be inspected weekly for inadequate water level. Inspections which give evidence of any particulation or discoloration suggest the water be replaced immediately.