



Recessed Showers, Eye Wash & Body Shower - mounting instruction

94G0110 - Rev. 10 May 2024



BROEN-LAB

Valve Cabinet - mounting instruction



17143009, 17143009-3289

17143909, 17143909-3288


Hole in wall, dimension

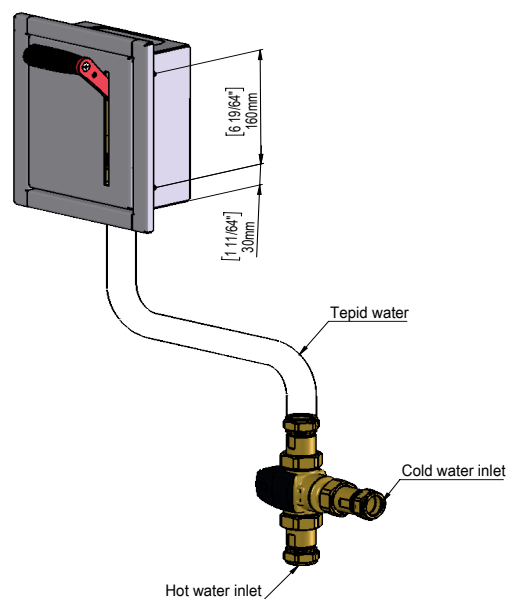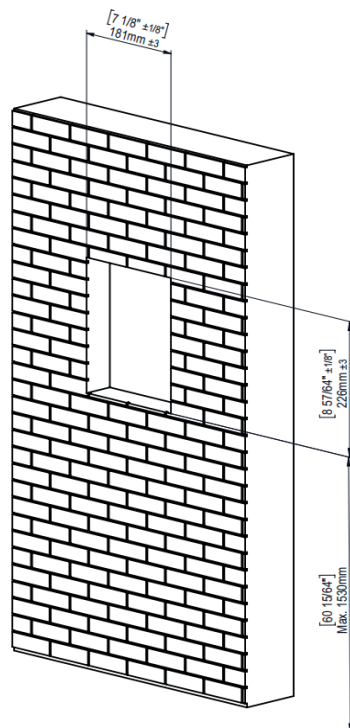

Cabinet, dimensions

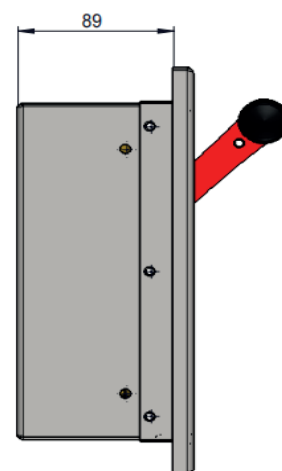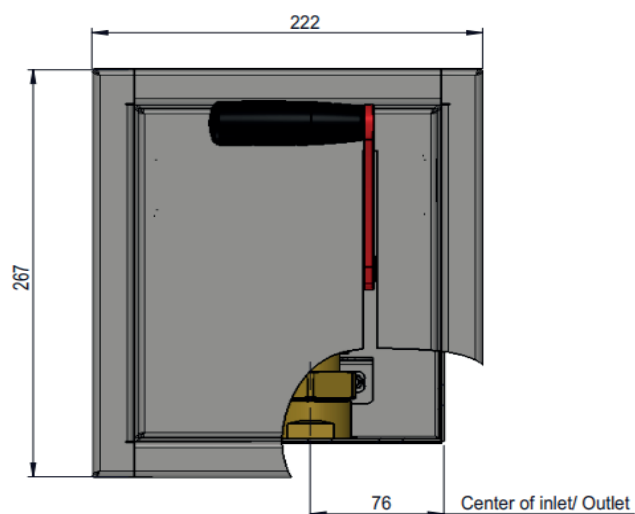

BROEN-LAB

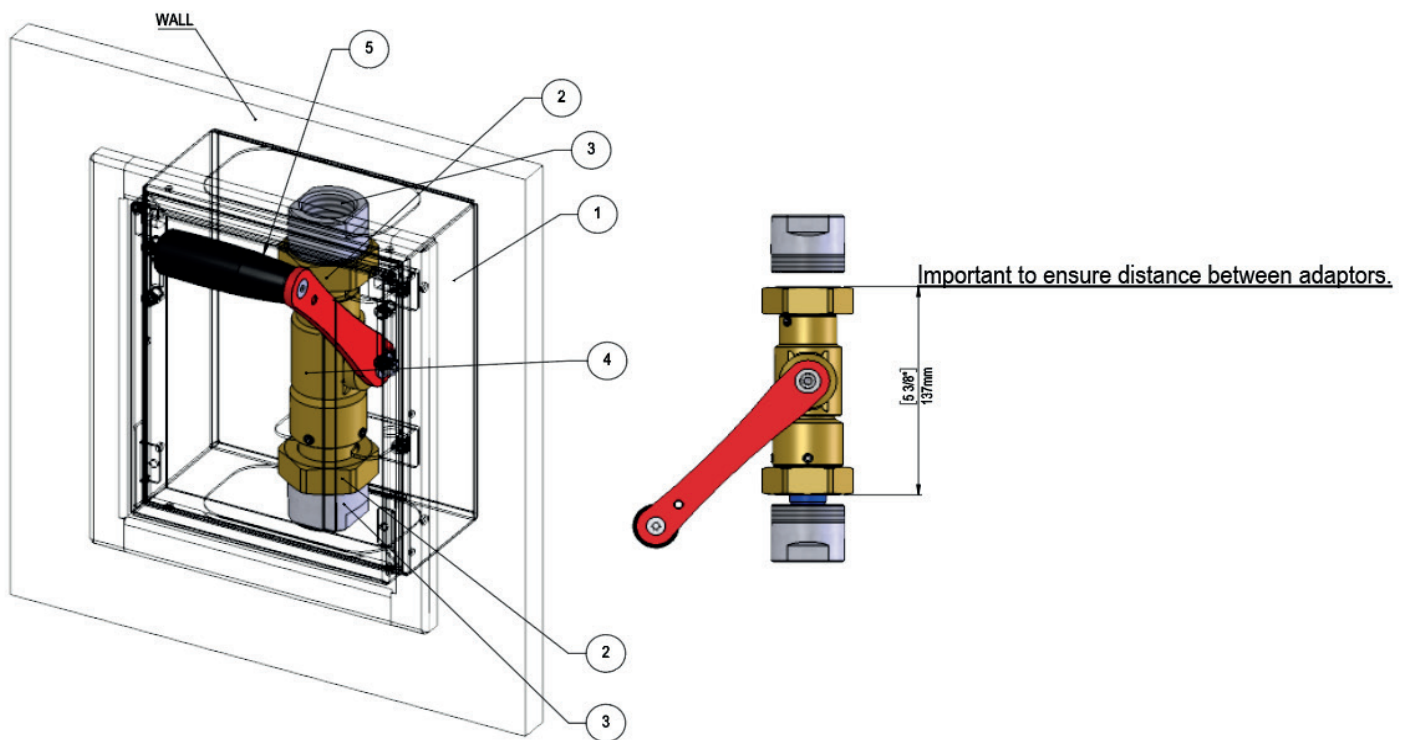

1. Attach the cabinet unit in the wall
2. Loosen the union nut for installing the pipe feed to the main pipe
3. Mount loose adaptors on the main pipe
4. Mount the body shower activator in bracket in cabinet and tighten union nut
5. Mount handle on shower activator arm



BROEN-LAB

Eye Wash - mounting instruction



17156009, 17156009-3289

17156909, 17156909-3289


Hole in wall, dimensions

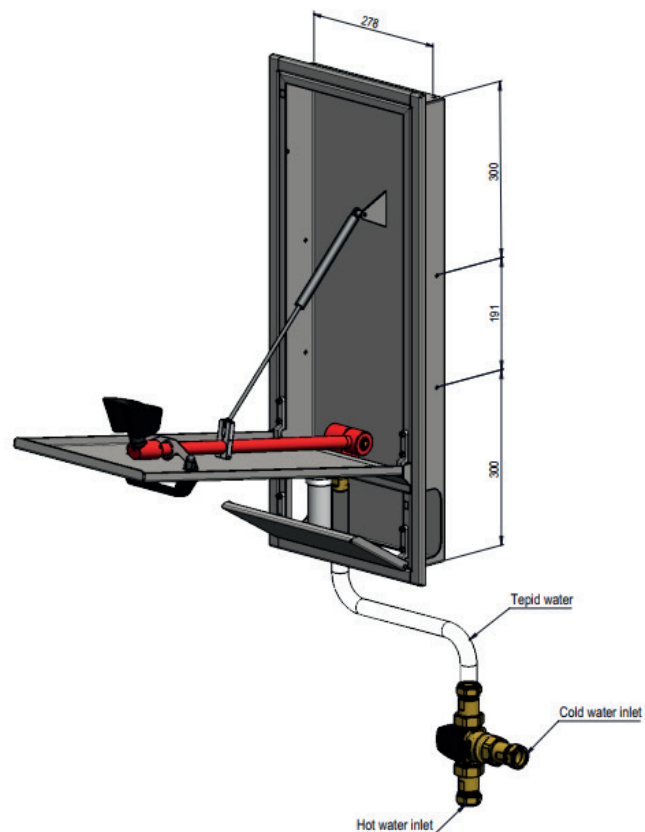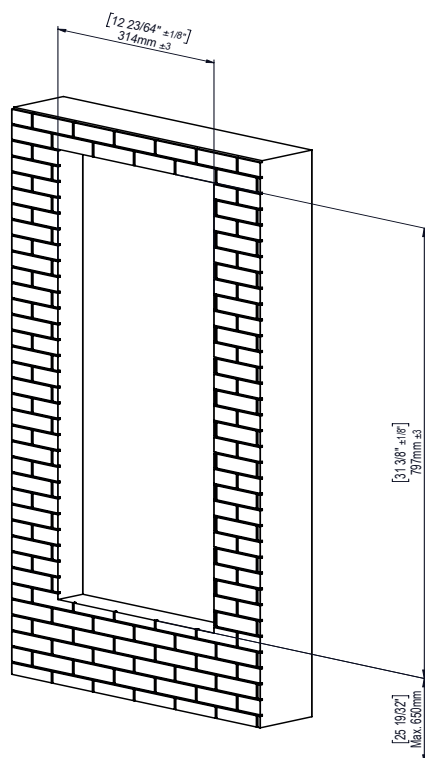

Cabinet, dimensions

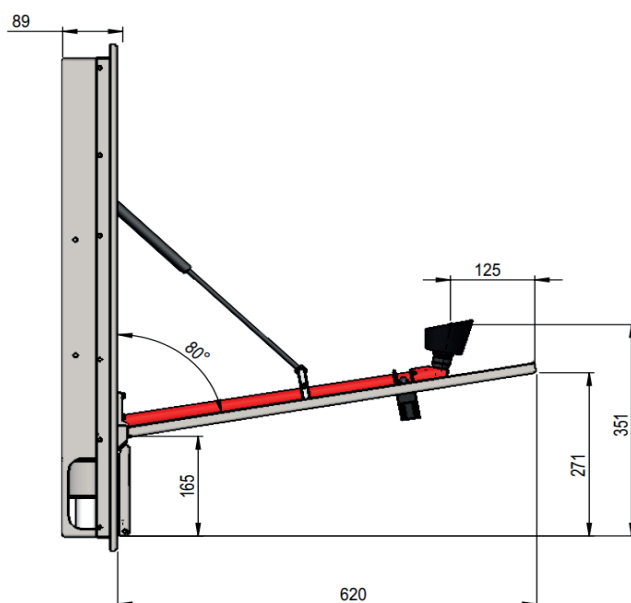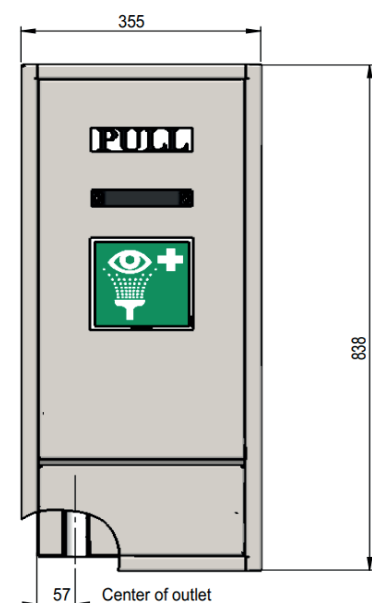

BROEN-LAB

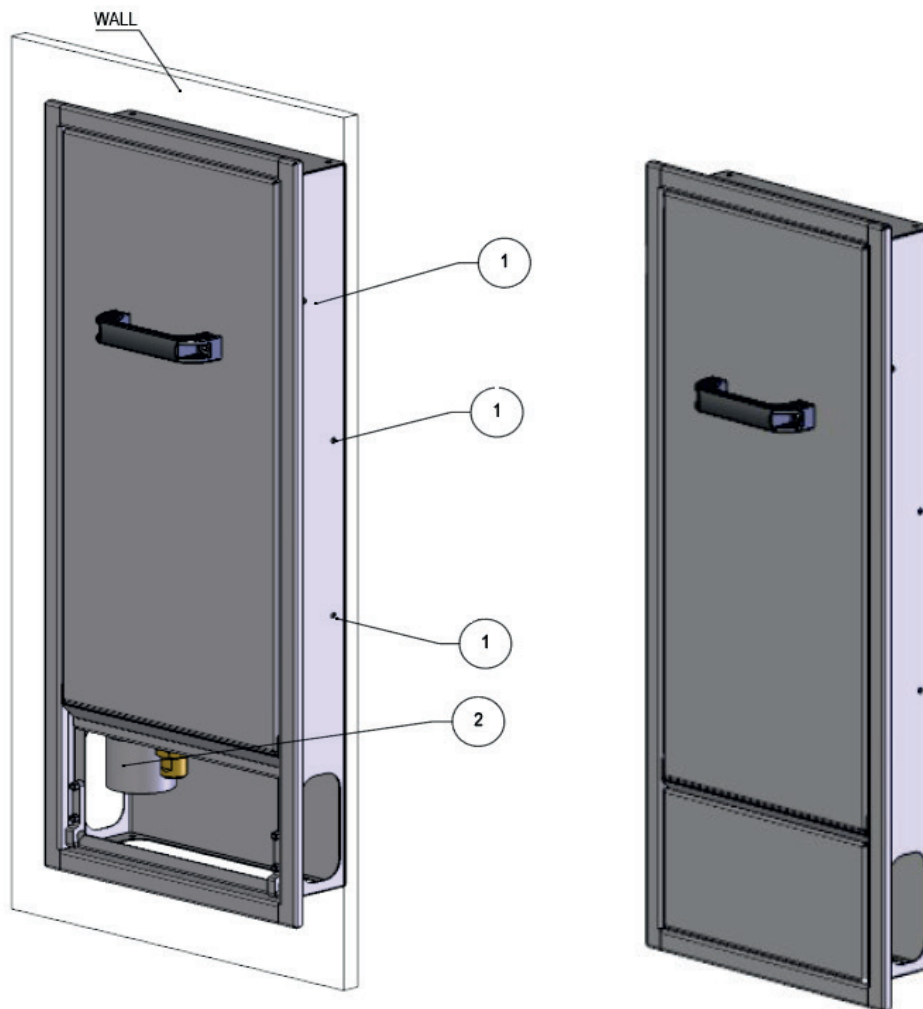

1. Attach the cabinet unit on both sides in the wall in the mounting holes
2. Connect the drain
3. Connect to the main pipe



BROEN-LAB

Eye Wash & Body Shower - mounting instruction



17157009, 17157009-3289

17157909, 17157909-3289

Hole in wall, dimension

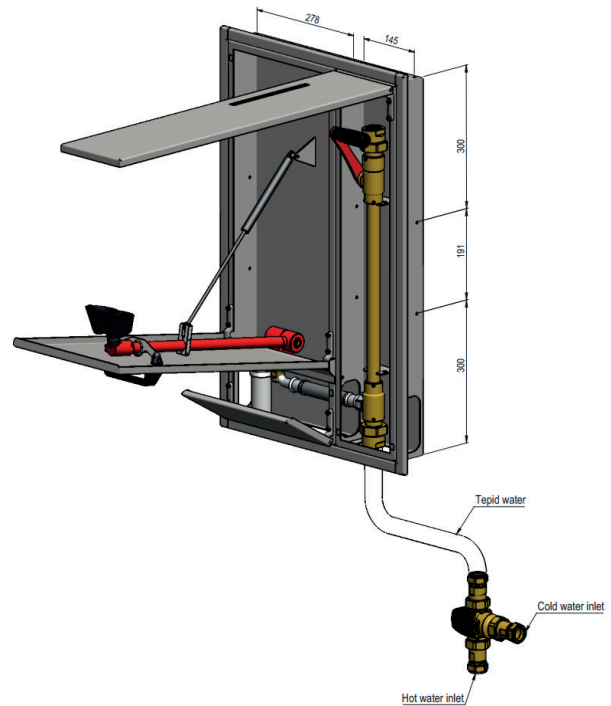
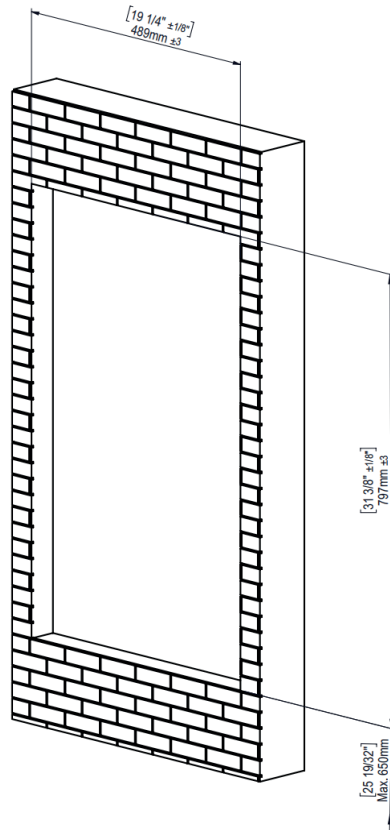


Cabinet, dimension

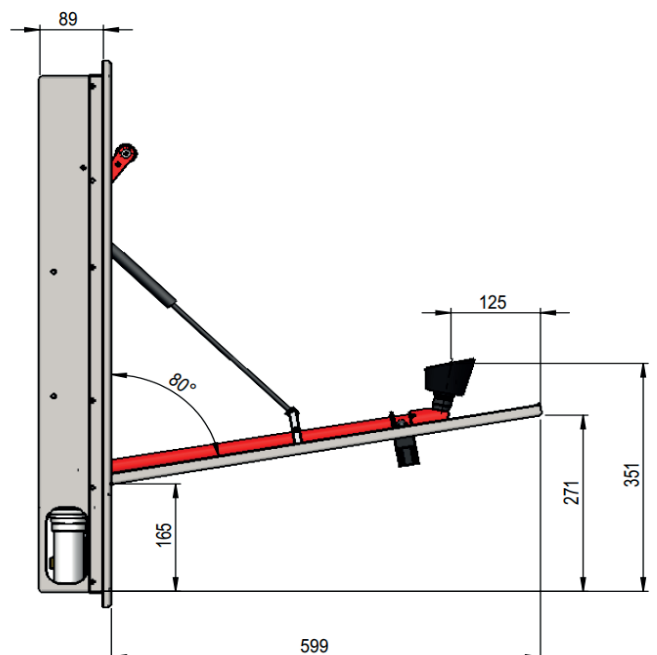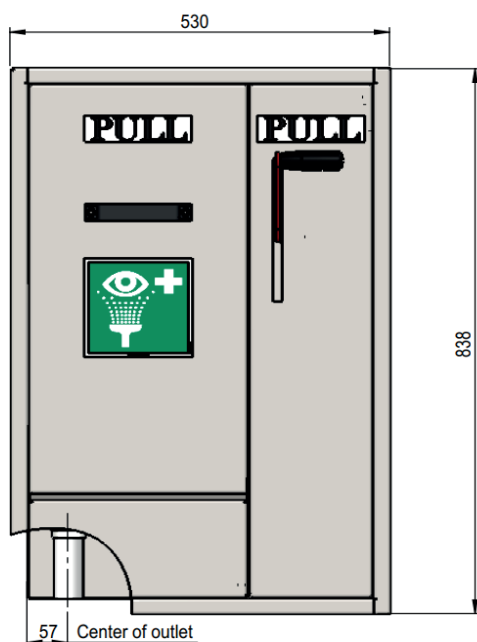

BROEN-LAB

BROEN-LAB

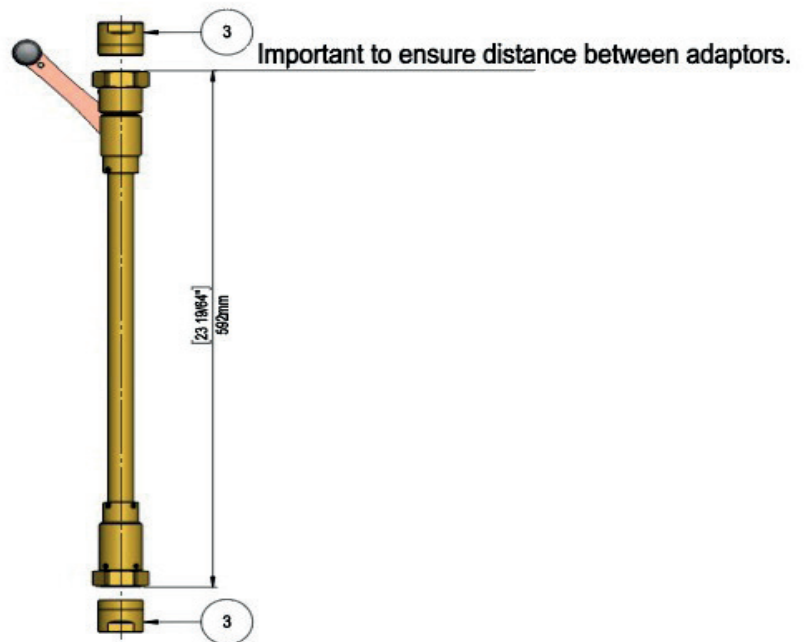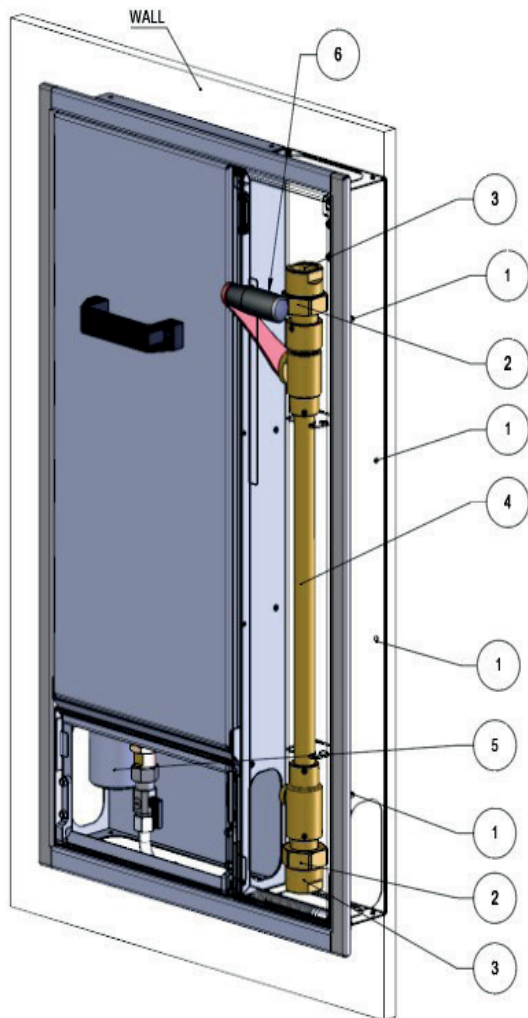

1. Attach the cabinet unit in the wall
2. Loosen the union nut for installing the pipe feed to the main pipe
3. Mount loose adaptors on main pipe
4. Mount the body shower activator in bracket in cabinet and tighten union nut
5. Connect the drain
6. Mount handle on shower activator arm

HEADQUARTER

BROEN-LAB A/S

Drejervaenget 2
5610 Assens
Denmark

Tel. +45 6376 6376

lab@broen-lab.com
broen-lab.com

INTERNATIONAL OFFICES

BROEN-LAB Ltd.

Tel. +44 121 522 4515

lab@broen-lab.com
broen-lab.com

BROEN-LAB Singapore Pte. Ltd.

25 International Business Park
#04-60A German Centre
Singapore 609916

Tel. +65 6298 0662

lab@broen-lab.com
broen-lab.com

BROEN-LAB GmbH

Tel. +45 6376 6376

lab@broen-lab.com
broen-lab.com

BROEN-LAB Sverige AB

Tel. +45 6376 6376

lab@broen-lab.com
broen-lab.com

BROEN-LAB Inc.

15 Constitution Drive
Suite 122
NH 03110
Bedford
USA

Tel. +1 603 310 5089

lab@broen-lab.com
broen-lab.com

BROEN-LAB develops, manufactures and sells Laboratory Fittings, Emergency Shower Systems and Eye Wash Systems into a broad spectrum of laboratories and industrial locations; sectors include Pharmaceutical, Food&Beverage and Academia. Our expertise and product quality ensure optimal solutions compliant to all relevant international norms setting new standards in risk mitigation in modern research and development facilities.

BROEN-LAB offers solutions that ensure the functionality, compliance, hygiene, durability and safety are of the highest priority. For more than 50 years our products have been integrated into a wide variety of workplaces, laboratories, hospitals and industrial locations, with features that are assessed and designed to be flexible, durable and compliant offering a broad range of options tailored to each location. This inherent design integrity provides confidence in the solution adopted wherever in the world these are applied.

BROEN-LAB is a collaboration partner in the design and layout of your next laboratory and it's integrated safety.



BROEN-LAB A/S ISO 9001 certification

In September 1991 BROEN-LAB A/S was certified according to ISO 9001 as one of the first Danish companies. The certification was carried out by Bureau Veritas, Denmark, for our Danish site in Assens.

The quality management system of BROEN-LAB A/S now complies with detailed specifications laid down by the internationally acknowledged EN ISO 9001:2015. This certification will further contribute to reputation for quality and reliability of BROEN-LAB A/S.

BROEN-LAB A/S ISO 14001:2015 certification

In December 2022 we got certified according to ISO 14001 in Environmental certification. The new certification was carried out by Bureau Veritas, Denmark, for Danish site in Assens. The Environment ISO14001 management system of BROEN-LAB A/S is included in internationally acknowledged EN ISO 9001:2015. Both certifications furthermore contribute to reputation for quality and reliability of BROEN-LAB A/S.